



## Details

### Ordering Information

For ordering please give **Article Number, Cap Cut, and Size**

(Example: CAP-96, C-Cut, 58):

- Article Number: **CAP-96**
- Cap Cut: **C-Cut** or **A-Cut**
- Size (given in cm head circumference):
  - Adult caps: **54, 56, 58, 60, 62, 64** (average male: 58, average female: 56)
  - Children caps: **52** (5-10 years), **54** (11-14 years)

The catalogue-number comprises the cap as described. For further information about accessories or consumables, please visit our website or contact our local distributor.

### Cap

White Subtemporal Cap with integrated chin belt

Sizes 52-64 made from High Precision Fabric

Options: C-Cut or A-Cut, Size. For further variations, please contact us.

The labelled holders are mounted according to the figure (1-96 transparent, Gnd black). The positions are chosen from 10%-System and from 5%-System<sup>1</sup>. Positions TP9/TP10 approximately meet left and right mastoids.

Each cap comes with 3 additional holders (apricot empty label) for use with double-sided adhesive rings to place electrodes on bare skin. The grip of the washer is better, when both, skin and holder, are degreased with alcohol prior to application.

### Accessories

The cap can be fixed either by the integrated chin belt alone, or additionally or alternatively by a chest belt (not included). Normally the decision between these 2 methods can be made according to what is more comfortable for the subject. Under difficult conditions the chest belt will hold the cap in place a little better.

### Service

In case of any further questions please contact our local distributor or [support@easycap.de](mailto:support@easycap.de). Please have the serial number of your cap at hand (syntax Axxxxx, located on the backside label of the cap).

---

<sup>1</sup> Electrode Nomenclature according to:

Oostenveld, R. & Praamstra, P. The five percent electrode system for high-resolution EEG and ERP measurements. *Clinical Neurophysiology* 2001; 112: 713-719