



EASYCAP

EEG Recording Caps and Related Products

EASYCAP GmbH

Am Anger 5

82237 Wörthsee

GERMANY

Tel +49 8153 88702-00

www.easycap.de

info@easycap.de

Standard 32Ch actiCAP Electrode Cap


Usable With and Without Dedicated Reference-Electrode

 Green Holders: Label 1-32, Hardware Channel 1 – 32

 Blue holder: Label & hard-wired Ref

 Black holder: Label & hard-wired Gnd

Electrode Names and Number Labels

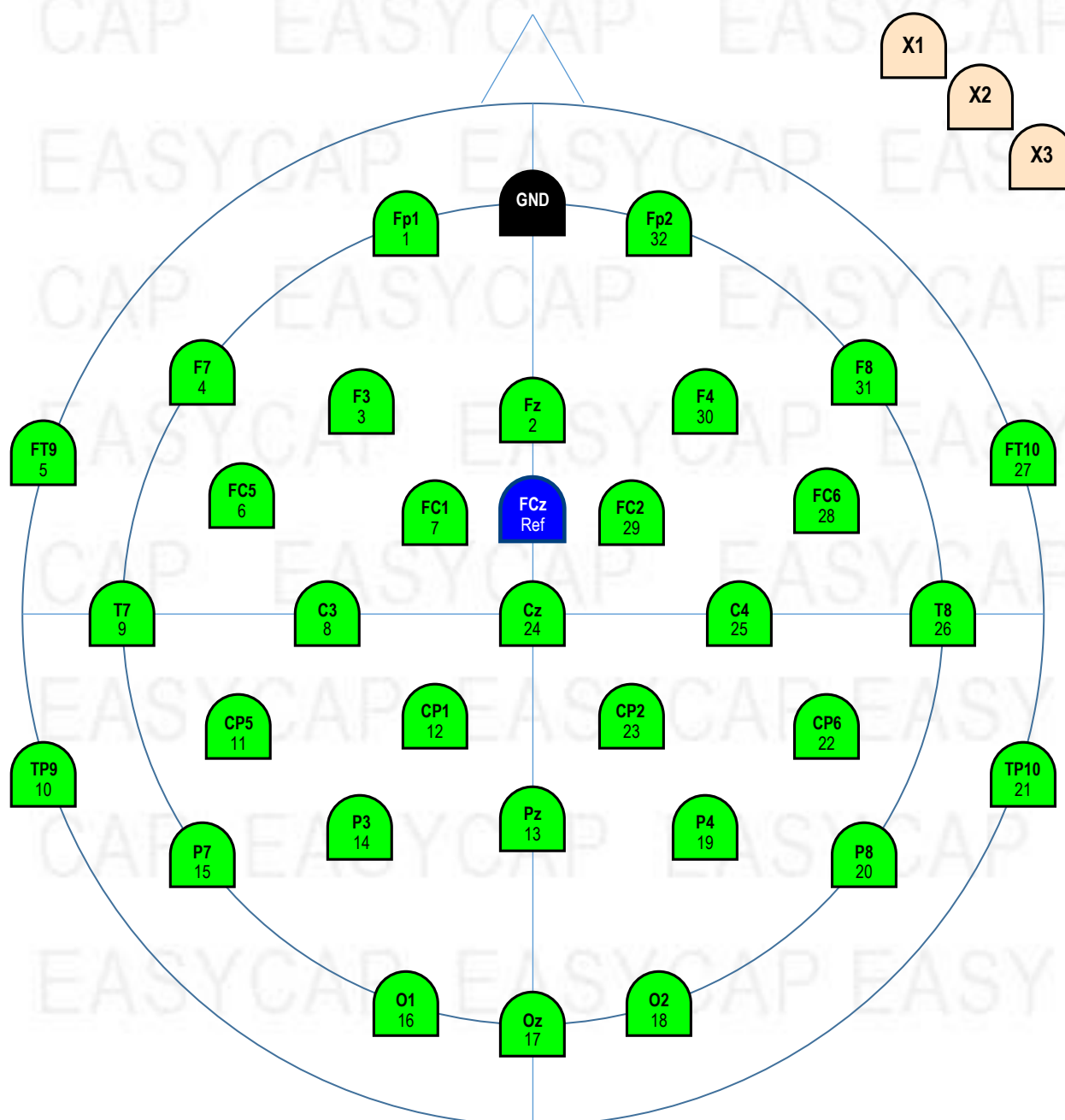

Details

Ordering Information

For ordering please give **Article Number, Cap Cut, and Size**
(Example: CMA-32, C-Cut, 58):

- Article Number: **CMA-32**
- Cap Cut: **C-Cut** or **A-Cut**
- Size (given in cm head circumference):
 - Adult caps: **52, 54, 56, 58, 60, 62, 64** (average male: 58, average female: 56)
 - Children caps: **50** (3-4 years), **52** (5-10 years), **54** (11-14 years)
 - Infant caps: **34, 36** (newborn), **38, 40** (3 months), **42, 44** (7 month), **46, 48** (2 years)

The catalogue-number comprises the cap as described. For further information about accessories or consumables please visit our website or contact our local distributor.

Cap

White SubInion Cap with integrated chin belt

Sizes 52 – 64 made from High Precision Fabric, Sizes 50 and smaller made from High Comfort Fabric

Options: C-Cut or A-Cut, Size. For further variations, please contact us.

The holders are mounted according to the figure. The positions include the 19 sites of the “International 10/20-System”. These core positions are extended by further positions of the 10%-System to increase the topographical resolution. Positions TP9/TP10 approximately meet left and right mastoids.

Each cap comes with 3 loose holders for use with double-sided adhesive rings to place electrodes on bare skin. The grip of the washer is better, when both, skin and holder, are degreased with alcohol prior to application.

Electrode Assignment

Plug the electrodes 1-32 into the green holders, which are also labelled 1-32. These are the amplifier channels 1-32.

If your amplifier requires a dedicated Ref-Electrode, plug Ref into the dark blue holder at FCz, otherwise leave the dark blue holder at FCz unused.

Supplementary Files

Each cap comes with a memory stick containing 2 bvef-files, one for use with REF, one for NO_REF. Use these bvef-files to assign the electrode names to channel numbers while creating the BrainVision Recorder Workspace. The bvef-file also contains the coordinates of the electrodes

Service

In case of any further questions

- Regarding consumables (e.g. electrolyte gel, syringes) and their usage, regarding cap, holders, layout, belts, please contact our local distributor or support@easycap.de. Please have the serial number of your cap at hand (syntax CMAxxxxx, located on the backside label of the cap).