



EASYCAP

EEG Recording Caps and Related Products

EASYCAP GmbH
Am Anger 5
82237 Würthsee
GERMANY


Tel +49 8153 88702-00
Fax +49 8153 88702-10
www.easycap.de
info@easycap.de

Standard 96Ch actiCAP Electrode Cap

Usable With and Without Dedicated Reference-Electrode

- Green Holders: Label 1-32, Hardware Channel 1 – 32
- Red Holders: Label 1-32, Hardware Channel 65 - 96
- Yellow Holders: Label 1-32, Hardware Channel 33 - 64
- Blue holder: For Ref or Ch64
- Black holder: Label & hard-wired Gnd

Electrode Names and Number Labels

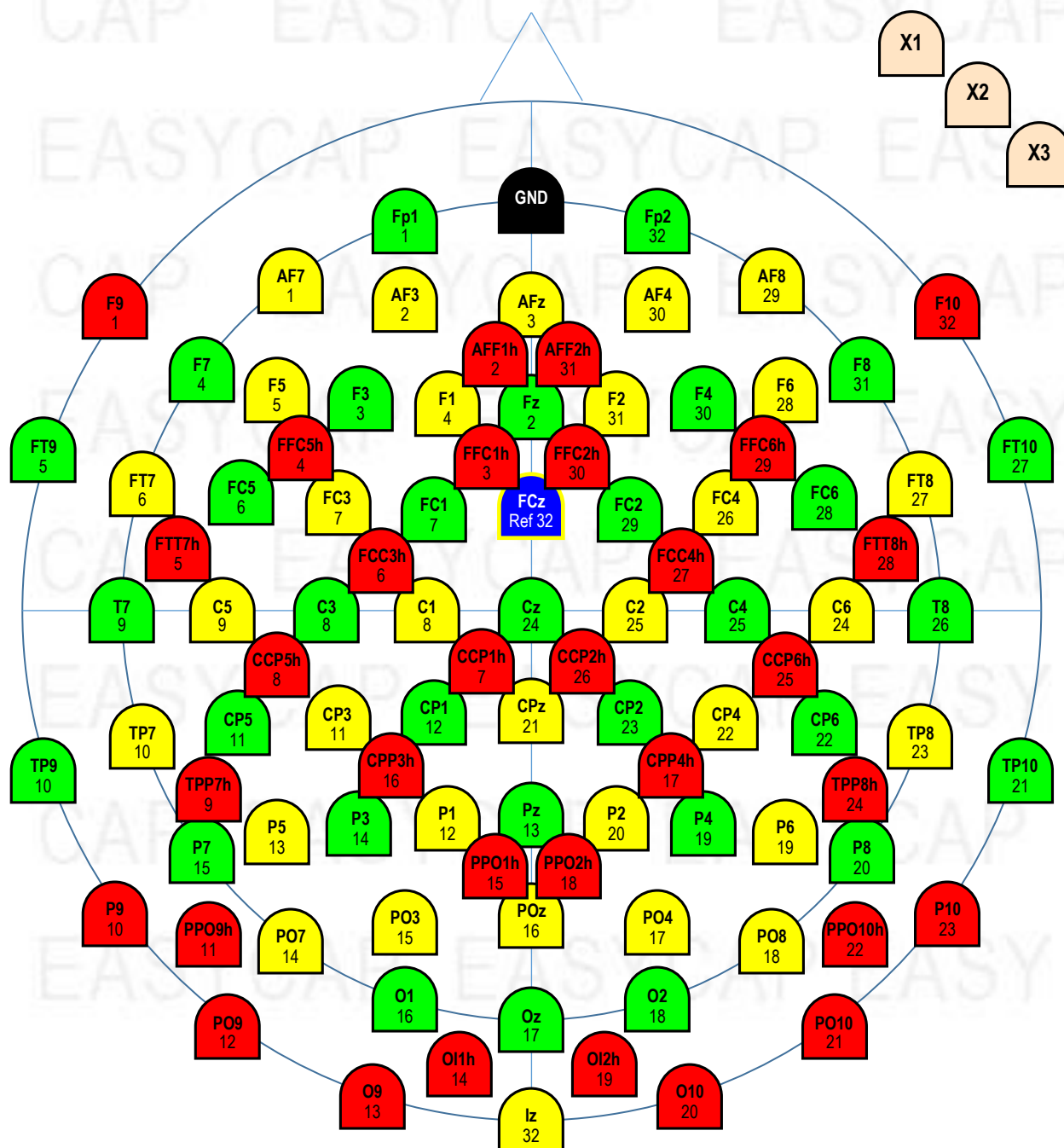

Details

Ordering Information

For ordering please give **Article Number, Cap Cut, and Size**
(Example: CMA-96, C-Cut, 58):

- Article Number: **CMA-96**
- Cap Cut: **C-Cut** or **A-Cut**
- Size (given in cm head circumference):
Adult caps: **52, 54, 56, 58, 60, 62, 64** (average male: 58, average female: 56)

The catalogue-number comprises the cap as described. For further information about accessories or consumables please visit our website or contact our local distributor.

Cap

White SubInion Cap with integrated chin belt, made from High Precision Fabric

Options: C-Cut or A-Cut, Size. For further variations, please contact us.

The holders are mounted according to the figure. The positions are chosen from 10%-System and from 5%-System¹. Positions TP9/TP10 approximately meet left and right mastoids.

Each cap comes with 3 loose holders for use with double-sided adhesive rings to place electrodes on bare skin. The grip of the washer is better, when both, skin and holder, are degreased with alcohol prior to application.

Electrode Assignment

The actiCAP electrodes come in bundles of 32, each labelled 1-32.

Plug the electrodes 1-32 of the first bundle into the green holders, which are also labelled 1-32. These are the amplifier channels 1-32.

Plug the second bundle into the yellow holders, which are channels 33-64. Red holders = Ch 65-96

If your amplifier requires a dedicated Ref-Electrode, plug Ref into the dark blue holder at FCz and plug Ch 64 (yellow 32) into Iz, otherwise plug Ch 64 (yellow 32) into the dark blue holder at FCz.

Supplementary Files

Each cap comes with a memory stick containing 2 bvef-files, one for use with REF, one for NO_REF. Use these bvef-files to assign the electrode names to channel numbers while creating the BrainVision Recorder Workspace. The bvef-file also contains the coordinates of the electrodes

Service

In case of any further questions

- Regarding consumables (e.g. electrolyte gel, syringes) and their usage, regarding cap, holders, layout, belts, please contact our local distributor or support@easycap.de. Please have the serial number of your cap at hand (syntax CMAxxxxx, located on the backside label of the cap).

WORKSHOP INTERNALS

¹ Electrode Nomenclature according to:

Oostenveld, R. & Praamstra, P. The five percent electrode system for high-resolution EEG and ERP measurements. Clinical Neurophysiology 2001; 112: 713-719