

Caution: To ensure correct and safe use of the products, make sure that you also read the Operating Instructions carefully prior to operation.



actiCAP CCCslim Setup




1 Installing the actiCAP software:



Plug in the USB stick and open the installation menu via double click on autorun.exe.

2 Important: Please make sure to respect the physical channel order when connecting electrode sets. Before connecting or disconnecting any electrode set (step 2 and 3) make sure that the actiCAP ControlBox is switched OFF.


Use the flat ribbon cable to connect the actiCAP ControlBox to the amplifier. Then connect each splitter box to the correspondent connector on the ControlBox.

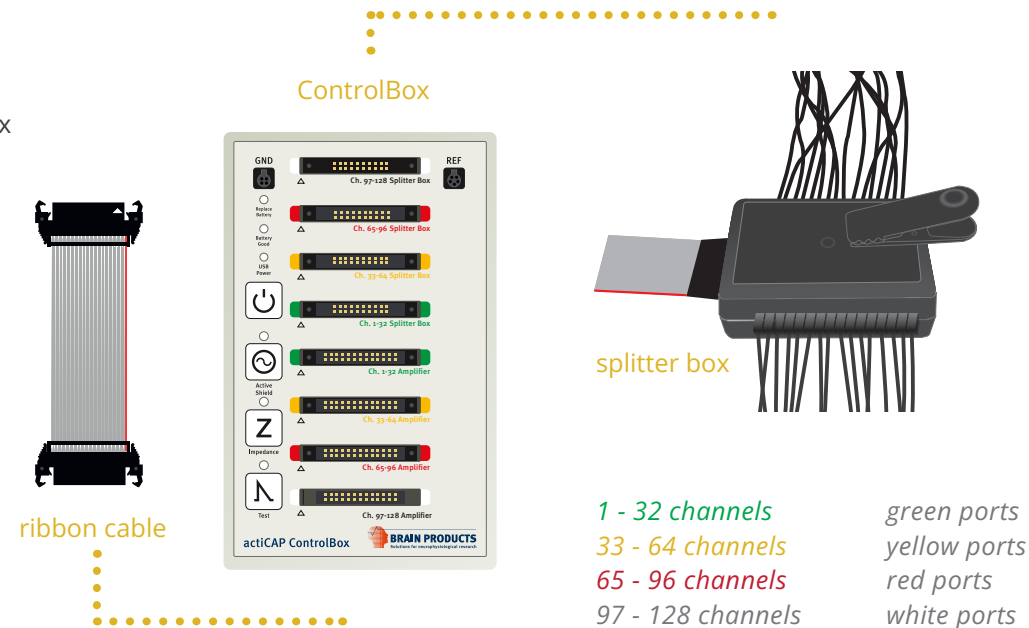


1 - 32 channels
33 - 64 channels
65 - 96 channels
97 - 128 channels

green ports
yellow ports
red ports
white ports

3 Connect the GND and REF electrode to the ControlBox

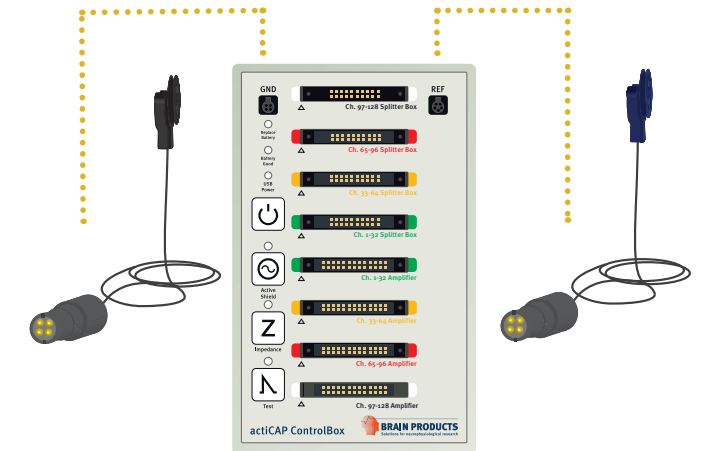


4 Installation of the actiCAP drivers

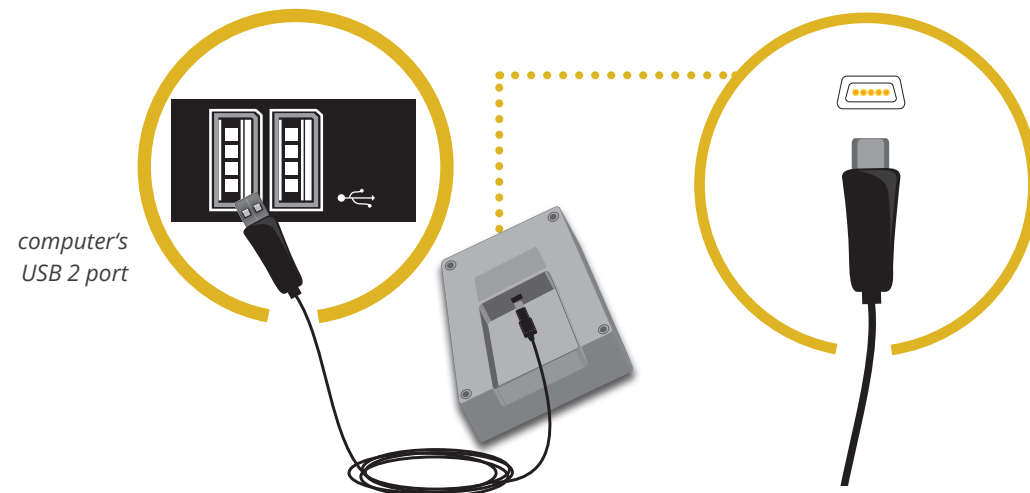

Remove the battery compartment cover on the rear side of the ControlBox, connect the actiCAP USB cable, and reattach the battery compartment cover. Then connect the actiCAP USB cable to your computer. Windows will install the driver automatically.

5

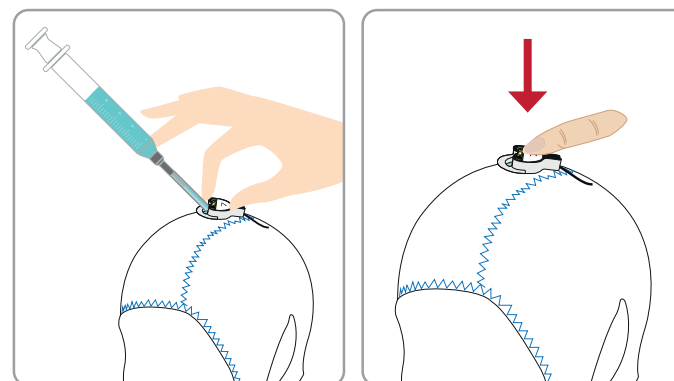



Prepare a 64 channel actiCAP slim or actiCAP snap electrode cap in approx. 10 minutes following these steps.

Start with the GND and REF electrodes.

- Slightly lift the electrode
- Inject the gel under the pellet
- Press down the electrode

Warnings:

- The splitter box must not come into contact with water.
- Do not pull the cables when removing the electrodes from the cap. Use both hands to remove the electrodes from the holder.

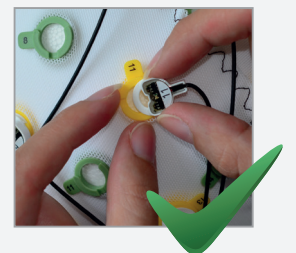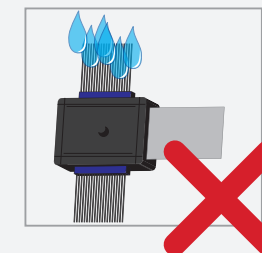

Setup Your actiCAP slim



Version 005
© Brain Products GmbH



Brain Products GmbH
Zeppelinstrasse 7
82205 Gilching
Germany
T +49 (0) 8105 733 84 0
F +49 (0) 8105 733 84 505
sales@brainproducts.com
www.brainproducts.com

Worldwide Caution:
Our products are scientific equipment for
INVESTIGATIONAL USE ONLY! Medical use
e.g. for diagnosis, treatment of disease or
other such purposes is strictly forbidden.